

List of Drug Concentrations for Various Cell lines to be used for Generating Stable Cell lines.

	<u>G418</u>	<u>Puro</u>	<u>Hygro</u>
HCT15	0.9mg/ml		400ug/ml
SAOS2	0.9mg/ml	1ug/ml	400ug/ml
U2OS	0.6mg/ml	1ug/ml	200ug/ml
IMEC	0.05mg/ml		25ug/ml

G418 – Media Tech (General Stores - Cat # 4361) Make up in 1x PBS. Store at –20°C.

Puro – Sigma (Cat #P7255) Store at –20°C.

Hygro – Boehringer Mannheim (Cat #843555) Store at 4°C.