

## Preparing Electroporation Cuvettes following use

- Rinse with distilled H<sub>2</sub>O (10X)
  - Dry over night
  - Incubate with 0.25N HCl for 30 min. (in glass beaker)
  - Wash with H<sub>2</sub>O, (5X)
  - Dry for several hours [upside down on Kimwipes (leaning against something so that they're at a slight angle)]
  - Incubate, with 95% ETOH in Tissue Culture Hood, 10 min.
  - Dry in Tissue Culture Hood
  - Put cap on
- That's it!